Test report

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Sample Information			
样品信息			
Sample Name	Nitrate	Sample properties	Solid
样品名称	硝酸根	样品性状	固体
Date of receipt	2025/11/11	During testing	2025/11/12 ~ 2025/11/14
收样日期		测试期间	
Test the ingredients and structural formula			
测试成分及结构式			
Nitrate (硝酸根)			
Experimental Requirements			
实验要求			
Analyze the nitrate solution in accordance with the customer's method, with the requirement that the retention time of nitrate remains stable at approximately 10 minutes. 按照客户方法分析硝酸盐溶液, 要求硝酸根出峰时间稳定在 10min 左右			
Reference Method			
参考方法			
Customer's method			
客户方法			
Reagent information			
试剂信息			
Reagent Name	Grade	Brand	Article Number
试剂名称	级别	品牌	货号

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1.2. Solution preparation (溶液配制)

1.2.1. Preparation of mobile phase (流动相配制):

Mobile phase: Weigh 4.28 g of potassium dihydrogen phosphate and dissolve it completely in 900 ml of purified water. Adjust the pH of the solution to 3.0 with phosphoric acid, mix well and filter. Combine the resulting solution with 100 ml of acetonitrile, mix thoroughly and sonicate to obtain the final solution.

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1.2.2. Blank solution、Diluent (空白溶液、稀释剂):

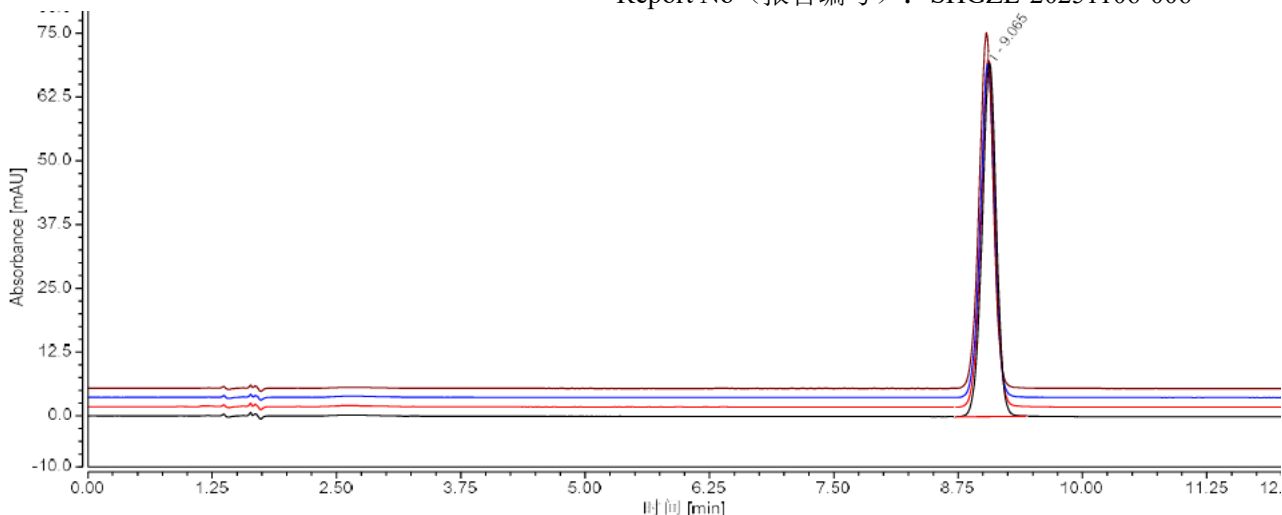
Methanol-Water (1:1, v/v) (等比例甲醇水)

1.2.3. Reference Solution: Weigh 3.6 mg of potassium nitrate reference substance into a 100 ml volumetric flask, dissolve it with solvent, dilute to the mark, and mix well to obtain the solution.

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
2. Chromatogram and data (谱图和数据)

(1) Analysis spectrum of Reference solution by column 1 (色谱柱 1 对照品溶液分析图谱):

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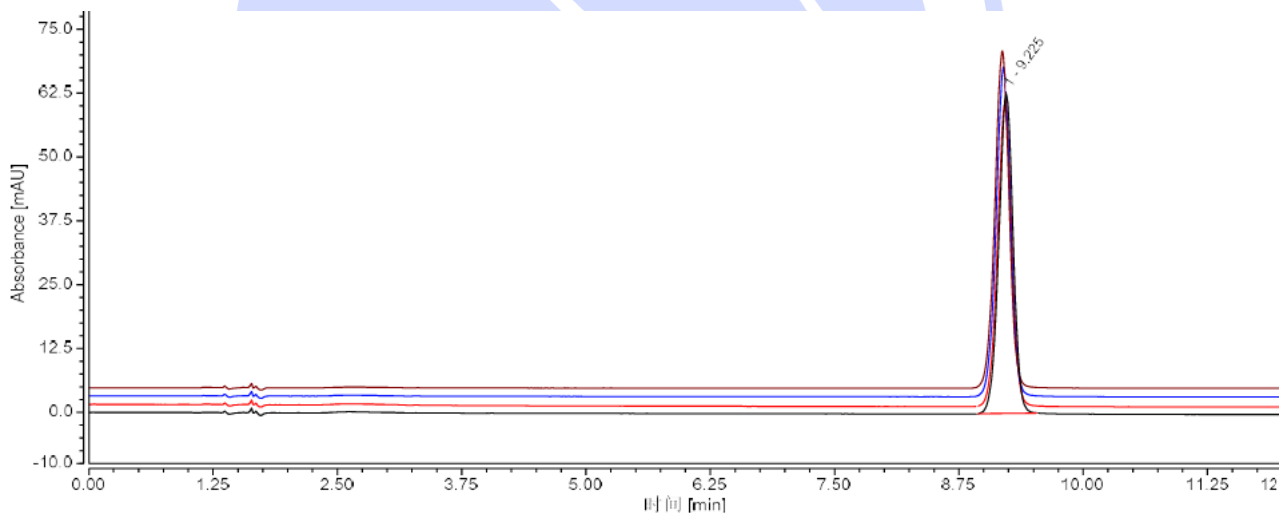(2) Analysis spectrum of Reference solution by coloum 2
(色谱柱 2 对照品溶液分析图谱):

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Web:www.welchmat.com

3. Conclusion (结论)

Using Welch Xtimate C18 (4.6×250mm, 5μm) analyzed reference solution under this chromatographic condition, the peak of the main component eluted at about 9min, which meets the customer's testing requirements.

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